

Rapid Isolation of RNA by Guanidinium Thiocyanate/Cesium Chloride Gradient Using Thermo Scientific Fiberlite Ultracentrifuge Rotors

Owen Mitch Griffith, PhD

KEY WORDS

- Cesium Chloride (CsCl) Gradient
- RNA Isolation
- Guanidinium Thiocyanate (GITC)

Introduction

The guanidinium thiocyanate method for isolation of ribonucleic acid (RNA) is used frequently when undegraded mRNA is desired⁽¹⁾. The RNA is pelleted through a high density cesium chloride (CsCl) solution to free it of DNA and polysaccharide contamination^(1,2). Customarily, this step is performed in a ultracentrifuge swinging bucket rotor for 12-22 hours. This report describes methods using Thermo Scientific Fiberlite carbon fiber fixed angle ultracentrifuge rotors to reduce run times to 4.5-5 hours allowing RNA isolation in a single day.

Procedures

The described RNA isolation protocol is derived from Sures and Crippa with minor modifications to prepare poly(A)+ RNA from cockroach (*Periplaneta americana*) heads^(3,4).

1. Harvest 25 g of tissue, freeze in liquid nitrogen (N₂) and grind to a powder under liquid N₂ with a mortar and pestle.
2. Resuspend powder in 150 mL of buffer containing 4.5 M guanidinium thiocyanate /50 mM EDTA (pH 8.0), 25 mM sodium citrate (pH 7.0), 0.1 M 2-mercaptoethanol, 2% lauroylsarcosine.
3. Homogenize suspension on a medium setting for 3-5 minutes at 4° C.
4. Remove insoluble material by centrifuging the suspension in a Fiberlite® F13-14x50cy carbon fiber rotor at 8,000 xg for 10 minutes at 4° C.
5. Add CsCl to the supernatant at a concentration of 0.2 g/ml.
6. Layer the sample solution over 10 mL of 5.7 M CsCl containing



Figure 1: Thermo Scientific Fiberlite F65L-6x13.5 rotor

50 mM EDTA (pH 7.3). If using a polyallomer tube, ensure the tube is completely full and the 5.7 M CsCl cushion should occupy one-third of the total tube volume. If necessary top off tube with 0.2 g/ml CsCl buffered solution.

7. Centrifuge sample solution in Fiberlite F50L-8x39 rotor at 50,000 rpm (~266,000 xg) for 5 hours at 20°C.

NOTE: Recommended conditions for other rotors are provided in Table 1.

8. After centrifugation, remove RNA pellet and dissolve in 2.0 mL of 10 mM Tris-HCl (pH 7.4), 2 mM EDTA.

9. To further purify, extract RNA 3 times with 2 volumes of phenol/chloroform (1:1) followed by 1 extraction with chloroform (Total RNA yield = 3.2 mg).

10. Isolate poly(A)+ RNA by 2 cycles of oligo(dT)-cellulose chromatography⁽⁵⁾ (Final poly(A)+ yield = 42 µg).

Conclusion

To confirm the integrity and biological activity of the RNA, the mRNA preparation can be translated *in vitro*. Using 200 ng of mRNA in a rabbit reticulocyte translation system, an 80 fold stimulation over background can be expected as determined by 35S-methionine incorporation into protein. The stimulation level will be comparable to that of the globin control. The translation products can be separated on a 12% polyacrylamide gel⁽⁶⁾. A wide range of protein sizes will be synthesized including some larger than 100kDa.

The mRNA can also be used in the construction of a cDNA library according to the first strand reaction conditions described by Gubler and Hoffman⁽⁷⁾. In previous studies, the yields were approximately 20% (µg cDNA/ µg mRNA) as judged by ³²P-labeled dCTP incorporation⁽⁷⁾. The cDNA products from this study had a wide size distribution including some

In addition to these offices, Thermo Fisher Scientific maintains a network of representative organizations throughout the world.



Figure 2: Thermo Scientific Fiberlite F50L-8x39 rotor

that appeared to be full-length copies of their corresponding mRNAs.

Using suggested protocols (Table 1), this method of RNA preparation in a Fiberlite carbon fiber fixed angle rotor, will result in a significant run time reduction. In order to maximize yield and stability use acceleration and deceleration rate of 8/8 (the second fastest setting) in Thermo Scientific Sorvall WX ultracentrifuges and the second fastest accel/dec rates in all other ultracentrifuges.

Rotor	Speed (rpm)	Time (h)
Fiberlite F50L-8x39	50,000	5.0
Fiberlite F65L-6x13.5	60,000	4.5

Table 1: Centrifugation conditions for RNA isolation

References

- Chirgwin, J.M., A.F. Przybyla, R.J. McDonald and W.J. Rutter, (1979). Isolation of biologically active ribonucleic acid from sources rich in ribonuclease. *Biochemistry* 18:5294-5299.
- Glisin, V., R. Crkvenjakov and C. Byus. (1974). Ribonucleic acid isolated by cesium chloride centrifugation. *Biochemistry* 13:2633-2637.
- Sures, I. and M. Crippa. (1984). Xenospin: the neurotensin-like octapeptide from *Xenopus* skin at the carboxyl terminus of its precursor. *Proc.Natl.Sci. USA.* 81:380-384.
- Turpen, T.H. and O.M. Griffith, (1986). Rapid Isolation of RNA

- by Guanidinium Thiocyanate/ Cesium Chloride Gradient Method. *BioTechniques* 4:1 13-14.
- Maniatis, T., E.F. Fritsch and J. Sambrook. (1982) *Molecular Cloning: A laboratory manual*. Cold Spring Harbor Lab., NY.
 - LeStourgeon, W.M. and A.L. Beyer. (1977) The rapid isolation, high-resolution electrophoretic characterization and purification of nuclear proteins, pp.387-406. In R. Stein, J. Stein and L. Kleinsmith (eds.), *Methods in Cell Biology*, Vol. 16. Academic Press, Inc., NY.
 - Gubler, U. and B.J. Hoffman. (1983). A simple and very efficient method for generating cDNA libraries. *Gene* 25:263-269.

North America:

USA / Canada
+1 866 984 3766
(866-9-THERMO)

Europe:

Austria
+43 1 801 40 0

Belgium
+32 2 482 30 30

France
+33 2 2803 2000

Germany national toll free
08001-536 376

Germany international
+49 6184 90 6940

Italy
+39 02 02 95059 434-254-375

Netherlands
+31 76 571 4440

Nordic/Baltic countries
+358 9 329 100

Russia / CIS
+7 (812) 703 42 15

Spain / Portugal
+34 93 223 09 18

Switzerland
+41 44 454 12 12

UK / Ireland
+44 870 609 9203

Asia:
China
+86 21 6865 4588 or
+86 10 8419 3588

India toll free
1800 22 8374

India
+91 22 6716 2200

Japan
+81 45 453 9220

Other Asian countries
+852 2885 4613

Countries not listed:
+49 6184 90 6940 or
+33 2 2803 2000

www.thermoscientific.com/centrifuge

© 2010 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.