

Accelerated Solvent Extraction (ASE) Techniques for In-Line Selective Removal of Interferences

Introduction

Interferences may be extracted along with desired analytes during an extraction process. These unwanted co-extractables may interfere with analyte detection or decrease instrument performance. Traditionally, chromatographic techniques such as gel-permeation chromatography (GPC) or a glass column packed with specific adsorbents are used to purify sample extracts prior to separation and analysis. Recent advances using Accelerated Solvent Extraction (ASE) systems, as described in several publications,^{1,3-10,12-17} include procedures for selective removal of interferences during sample extraction, thus combining extraction and purification into a single step.

This application note summarizes seven ASE procedures developed to remove co-extractable material from various matrices: procedures to selectively extract polar compounds from lipid-rich samples and to fractionate lipids from biological samples.

This note is intended to serve as a guide to develop ASE methods. For more information, please refer to the original publication cited with each method described below, or contact us.

Selective Extraction of Nonpolar Compounds

In an effort to eliminate post-extraction cleanup steps, we and others have researched the addition of various adsorbents to the extraction cell. For many sample types, this approach has proven successful in producing clean extracts that are ready for direct analysis. For example, nonpolar lipids are often co-extracted from fish tissue. Adding alumina (aluminum oxide, Al_2O_3 , acidic, activated by placing in a drying oven at 350 °C for 15 h) to the extraction cell before adding the sample or sample mixture has been shown to prevent the extraction of unwanted lipids. Mixing the sample with C18 resin (1:2) has been shown to retain organic contaminants. (C18 bonded silica, 35–70 μm diameter, and porosity of 60 Å, from Alltech has been used, but similar materials from other vendors can be used.)

When in-cell cleanup is performed during ASE, solvent choice impacts the retention of unwanted components. For example, a mixture of hexane/acetone (1:1) is a common solvent for extracting organochlorine pesticides from animal tissues. After extraction, a cleanup step is usually required to remove co-extracted lipids. Adding alumina to an ASE cell and extracting with hexane can prevent the extraction of interferences. However, if hexane/acetone (1:1) is used as the extraction solvent, almost no lipid material will be retained on the alumina. Table 1 lists two types of fat retainers and the ratio of each required to retain fat when using nonpolar solvents. Table 2 lists common adsorbents that are used for selective extraction of compounds using ASE systems.

Selected Techniques for In-Line Cleanup of Various Nonpolar Compounds

Selective Extraction of PCBs from Fish Meal Samples Using Acid-Impregnated Silica Gel^{5,6} Adsorbents and Dispersants

Silica gel
Thermo Scientific Dionex ASE Prep DE (P/N 062819)
Fisher Scientific sodium sulfate*
Fisher Scientific Ottawa sand
* Sodium sulfate can be used only when using hexane or heptane as an extraction solvent. If using any other solvent, it is important to use Dionex ASE™ Prep DE.

Table 1. Ratio of Fat Retainer per Gram of Lipid Using Nonpolar Solvents

Fat Retainer	Ratio of Fat Retainer/Lipid (g/g)
Alumina	20–25
Acid-impregnated silica gel	45.5

The nonpolar solvents that give the best results are hexane and heptane. Increased extraction temperatures and increased solvent polarity will decrease the fat retainer's effectiveness.

Table 2. List of Adsorbents and Uses

Adsorbent*	Uses
Carbon	Removes organics and nonpolar compounds (used in dioxin analysis) ¹
Copper	Removes sulfur ²
Ion-exchange resins**	Removes organics, ionic interferences for IC and IC/MS analysis ³
CS18 resin	Removes organics, polar compounds, lipids, colors ⁴
Acid-impregnated silica gel**	Removes lipids ^{5,6}
Alumina**	Removes nonpolar lipids, colors ^{7,8,9}
Florisil® (U.S. Silica Co., Frederick, MD, USA)	Removes nonpolar lipids ^{10,11}
Silica gel	Removes nonpolar lipids ¹²

**Note: We do not recommend using sodium sulfate and magnesium sulfate as in-cell adsorbent agents with polar solvents because of potential clogging problems. Use of sodium or magnesium sulfate can also contribute to the failure of key ASE instrument components such as the static valve.

**These selected techniques are detailed below.

Preparation of 40% Sulfuric Acid-Impregnated Silica Gel

Heat 600 g of silica gel at 200 °C for at least 12 h. Allow to cool. Add 400 g of sulfuric acid to the cooled silica gel. Mix together in a tumbling mixer for 4 h.

ASE Extraction Conditions

Extraction Solvent:	n-Hexane
Pressure:	1500 psi
Temperature:	100 °C
Static Time:	5 min
Static Cycles:	2
Flush:	60%
Purge:	90 s
Cell Size:	33 mL

Extraction Cell Preparation

Prior to extraction, prepare a 33-mL cell as shown in Figure 1. Begin by inserting two cellulose filters at the outlet end of the ASE cell. Add a layer of untreated silica gel then a layer of the acid-impregnated silica gel. Place a cellulose filter on top of the acid-impregnated silica gel then add the sample, as prepared below. Fill any remaining cell volume with Ottawa sand. (Placing a cellulose filter on top of the Ottawa sand is optional.)

Sample Preparation

Grind 2 g of fish meal sample with 2 g of sodium sulfate and 2 g Ottawa sand using a mortar and pestle. Add this sample mixture to the cell as shown in Figure 1. Place the extraction cell, containing the sample, on the ASE and extract with the method conditions listed above. Concentrate the extracts to approximately 1 mL then dilute to a final volume of 2 mL with hexane. Analyze using EPA Method 8082.

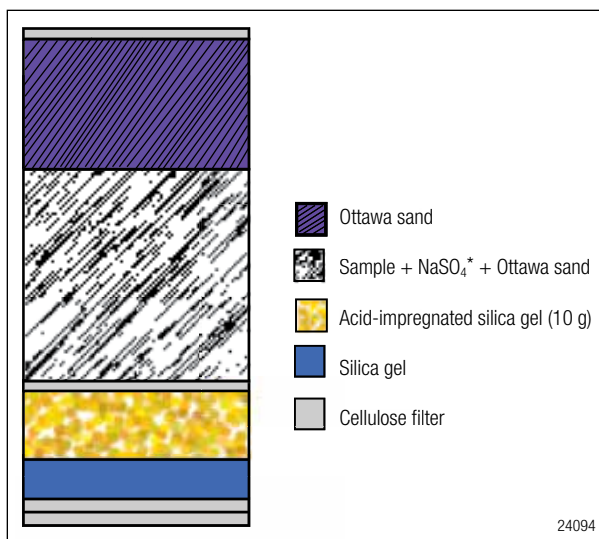


Figure 1. Preparation of ASE cell for the selective extraction of PCBs from fish meal.

Extraction Cell Preparation

Prior to extraction, prepare the 33-mL cell by placing one cellulose filter at the outlet. Place the activated alumina on top of the filter, as shown in Figure 2.

Sample Preparation

Grind 3 g of fish tissue with 15 g of sodium sulfate in a mortar and pestle. Transfer this mixture to the 33-mL extraction cell as shown in Figure 2. Rinse the mortar and pestle with 2 mL of hexane and add to the extraction cell. Place the extraction cell containing the sample on the ASE and extract with the method conditions listed above. Analyze extracts using EPA Method 8082.

The amount of alumina required to retain the fat in the sample depends on the fat content of the fish tissue and the mass of the tissue sample extracted (Table 1).

Selective Extraction of PCBs from Fish Tissue Using Alumina⁷

Adsorbents and Dispersants

Dionex ASE Prep DE (P/N 062819)

Fisher Scientific sodium sulfate*

Alumina 60–325 mesh, Brockman activity I, Fisher Scientific acidic

*Sodium sulfate can be used only when using hexane or heptane as an extraction solvent. If using any other solvent, it is important to use Dionex ASE Prep DE.

Preparation of Activated Alumina

Heat the alumina at 350 °C for 15 h.

ASE Extraction Conditions

Extraction Solvent:	n-Hexane
Pressure:	1500 psi
Temperature:	100 °C
Static Time:	5 min
Static Cycles:	2
Flush:	60%
Purge:	120 s
Cell Size:	33 mL

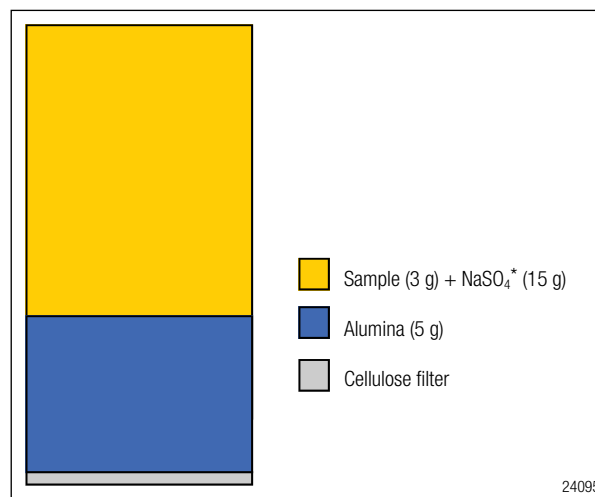


Figure 2. Preparation of ASE cell for the selective extraction of PCBs from fish tissue using alumina.

Selective Extraction of Polar Compounds from Polar Matrices

Selective Extraction of Perchlorate from Vegetation Using Ion-Exchange Resins^a

Ion-Exchange Resins and Adsorbents

Thermo Scientific Dionex OnGuard II Sample Pretreatment Cartridges

Ag (P/N 057089)

Ba (P/N 057093)

H (P/N 057085)

RP (P/N 057083)

Dionex ASE Prep DE (P/N 062819)

Fisher Scientific basic alumina

ASE Extraction Conditions

Extraction Solvent:	HPLC-grade water
Pressure:	1500 psi
Temperature:	80 °C
Static Time:	5 min
Static Cycles:	3
Flush:	30%
Purge:	120 to 240 s
Cell Sizes:	33 mL or 100 mL

Extraction Cell Preparation

Prior to extraction, prepare the 100-mL cell by inserting two glass-fiber filters in the outlet end of the ASE cell. Cut the end off each Dionex OnGuard™ resin cartridge and store the resins in separate beakers until needed. Add the OnGuard resins, the alumina, and the glass-fiber filters as shown in Figure 3. Fill any remaining cell volume with Dionex ASE Prep DE. If using 33-mL cells, prepare in the same manner with proportionally less of each adsorbent. To ensure clean adsorbents and resins, extract each prepared cell with the ASE method conditions listed above before adding the sample.

Sample Preparation

Remove 10 g of “clean” Dionex ASE Prep DE from the extraction cell and grind together with 5 g of vegetation sample. Add the sample mixture back to the extraction cell. Place the extraction cell containing the sample on the ASE system and extract with the method conditions listed above. Adjust the final volume of each extract to either 40 mL (if a 33-mL ASE cell was used) or 100 mL (if a 100-mL cell was used).

Selective Extraction of Corticosteroids from Bovine Liver¹³

This method describes a fractionation extraction to remove nonpolar triglycerides (defatting step) followed by a second extraction of the sample to recover analytes of interest.

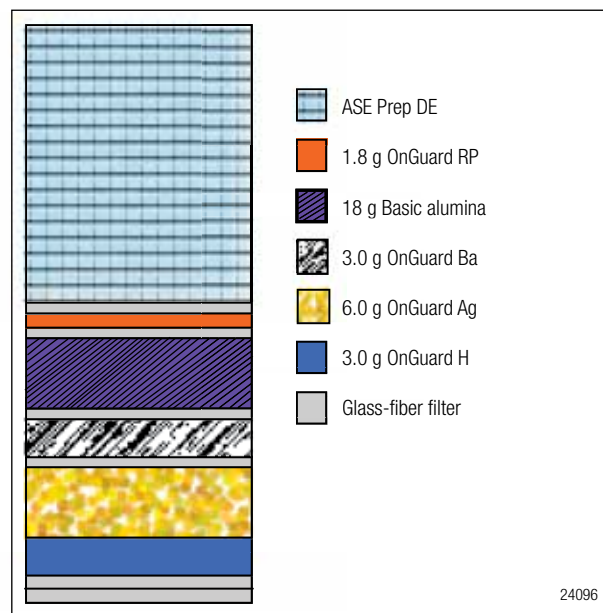


Figure 3. Preparation of ASE cell for the selective extraction of perchlorate from vegetation.

ASE Extraction Conditions

	Defatting	Extraction
Solvent:	n-Hexane	Hexane/ethyl acetate (1:1, v/v)
Temperature:	60 °C	50 °C
Pressure:	1500 psi	1500 psi
Static Time:	5 min	5 min
Flush:	100%	60%
Purge:	100 s	100 s
Static Cycles:	3	1
Cell Size:	22 mL	22 mL

Sample Preparation

Mix 5 g of liver tissue sample with 7.5 g of Dionex ASE Prep DE and load into a 22-mL extraction cell containing a cellulose filter at the outlet end. First extract with the defatting conditions listed above and re-extract the sample using the analyte extraction conditions. Evaporate second extract to dryness and dissolve in 4 mL of absolute ethanol. Centrifuge at 4000 rpm. Remove the ethanol layer and evaporate to dryness. Dissolve residue in 500 µL of methanol. Analyze by LC/MS.

Selective Extraction of Sulfonamide from Animal Tissue^{4,14}

This method describes the use of C18 resin to remove most of the nonpolar lipids from animal tissue samples during extraction. The remaining co-extracted polar lipids are precipitated from the extract by cooling and separated by centrifugation. The resulting supernatant is analyzed using HPLC/MS/MS.

ASE Extraction Conditions

Extraction Solvent:	HPLC-grade water
Pressure:	1500 psi
Temperature:	160 °C
Static Time:	5 min
Static Cycles:	1
Flush:	60%
Purge:	60 s
Cell Size:	11 mL

Sample Preparation

Homogenize the meat samples using any standard tissue homogenizer. This should be done with HPLC-grade water added to the sample. Set the homogenizer to 5,000 rpm. Once the homogenizer reaches 5,000 rpm, increase speed to 25,000 rpm and hold for 15 min. Evaporate excess water. Weigh out approximately 2 g of homogenized tissue and mix with 4 g of C18 until the entire mixture is of a uniform consistency. Transfer this mixture to an 11-mL ASE cell containing a glass-fiber filter.

Place the cell onto the Thermo Scientific Dionex ASE 200 instrument. Program the Dionex ASE instrument with the conditions above and start the extraction. After the extraction, place the sample vials in a freezer (–18 °C) for one hour to precipitate the lipids.¹⁴ Remove the extracts from the freezer and centrifuge for 5 min at 10,000 rpm. Analyze the supernatant using HPLC/MS/MS as described in reference 4.

Extraction of Acrylamide from Cocoa and Coffee Samples Using In-Cell Cleanup¹⁵

The addition of Florisil to the extraction cell greatly simplifies the analysis of acrylamide from complex food samples such as cocoa and coffee. This fact is illustrated in Figure 4. Coffee samples, each with different amounts of Florisil added to the ASE cell, were extracted and loaded into Isolute[®] (Biotage AB, Uppsala, Sweden) Multimode SPE cartridges. Figure 4 clearly shows a decrease in the amount of visible co-extractable material as the amount of Florisil in the ASE cell increased. The authors report that 6 g of Florisil is sufficient to produce a clear extract. For complete details of the ASE conditions and sample preparation see Application Note 358.¹⁵

Lipid Fractionation of Biological Samples^{16,17}

The following ASE method was developed to fractionate nonpolar neutral lipids and polar phospholipids from biological samples. The fractionation takes place in a single ASE cell containing a specific adsorbent that retains polar phospholipids while nonpolar lipids are extracted. Once the nonpolar lipids are extracted, the sample is re-extracted with a more polar solvent, which removes the polar phospholipids. Each extract is directed into a separate collection vial, allowing fractionation of the two classes of lipids. This fractionation is feasible due to the flow-through design of the ASE. The solvent controller allows the sequential use of different extraction solvents. For example, a single ASE schedule can be used to automate an extraction with hexane/acetone (9:1) at 50 °C to remove triglycerides, followed by an extraction with chloroform/methanol (1:4) to isolate polar lipids such as phospholipids or hydroxyl-fatty acids. Additional details about this method can be found in references 16 and 17.

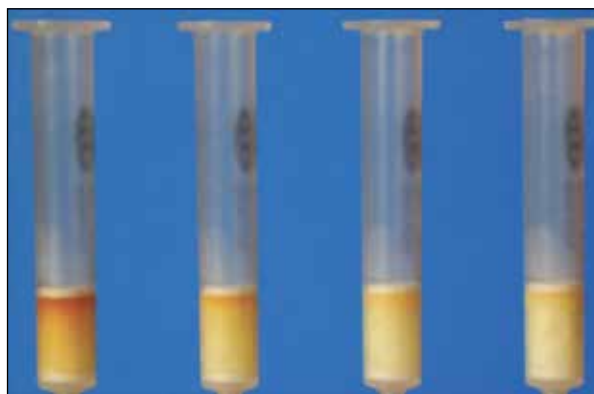


Figure 4. Residual co-extracts after the ASE extraction of a coffee sample trapped on an Isolute Multimode SPE cartridge, the ASE cells containing from 0 to 6 g Florisil. From left to right: 1) no Florisil in the ASE cell, 2) 2 g Florisil in the ASE cell, 3) 4 g Florisil in the ASE cell, and 4) 6 g Florisil in the ASE cell.

Conclusion

The flow-through design of ASE allows the system to perform extractions in combination with in-line cleanup. Eliminating post-extraction cleanup steps reduces solvent usage and waste, and saves the operator time and labor. This makes the ASE system a powerful tool for total sample preparation.

References

- Nording, M.; Sporning, S.; Wiberg, K.; Bjorklund, E.; Haglund, P. Monitoring Dioxins in Food and Feedstuffs Using Accelerated Solvent Extraction with a Novel Integrated Carbon Fractionation Cell in Combination with CAFLUX Bioassay. *Anal. Bioanal. Chem.* **2005**, *381*, 1472-1475.
- U.S. EPA Method 3660B, Sulfur Cleanup. U.S. Environmental Protection Agency, Cincinnati, OH, 1996.
- Dionex Corporation (now part of Thermo Fisher Scientific). Determination of Perchlorate in Vegetation Samples Using Accelerated Solvent Extraction (ASE) and Ion Chromatography. Application Note 356; LPN 1807. Sunnyvale, CA, 2006.
- Dionex Corporation (now part of Thermo Fisher Scientific). Rapid Determination of Sulfonamide Residues in Animal Tissue and Infant Food Containing Animal Products Using Accelerated Solvent Extraction (ASE). Application Note 353; LPN 1708. Sunnyvale, CA, 2005.
- Bjorklund, E.; Muller, A.; von Holst, C. Comparison of Fat Retainers in Accelerated Solvent Extraction for the Selective Extraction of PCBs from Fat-Containing Samples. *Anal. Chem.* **2001**, *73*, 4050-4053.
- Sporring, S.; Bjorklund, E. Selective Pressurized Liquid Extraction of Polychlorinated Biphenyls from Fat-Containing Food and Feed Samples Influence of Cell Dimensions, Solvent Type, Temperature and Flush Volume. *J. Chromatog., A.* **2004**, *1040*, 155-161.
- Ezzell, J.; Richter, B.; Francis, E. Selective Extraction of Polychlorinated Biphenyls from Fish Tissue Using Accelerated Solvent Extraction. *American Environmental Laboratory.* **1996**, *12*, 12-13.
- Dionex Corporation (now part of Thermo Fisher Scientific). Selective Extraction of PCBs from Fish Tissue Using Accelerated Solvent Extraction (ASE). Application Note 322; LPN 0764. Sunnyvale, CA, 1996.
- Dionex Corporation (now part of Thermo Fisher Scientific). Determination of PCBs in Large-Volume Fish Tissue Samples Using Accelerated Solvent Extraction (ASE). Application Note 342; LPN 1204. Sunnyvale, CA, 2000.
- Gomez-Ariza, J.L.; Bujalance, M.; Giraldez, I.; Velasco, A.; Morales, E. Determination of Polychlorinated Biphenyls in Biota Samples Using Simultaneous Pressurized Liquid Extraction and Purification. *J. Chromatog., A.* **2002**, *946*, 209-219.
- U.S. EPA Method 3620C, Florosil Cleanup. U.S. Environmental Protection Agency, Cincinnati, OH, 2000.
- Gomez-Ariza, J.L.; Garcia-Barrera, T.; Lorenzo, F.; Gustavo Gonzales, A. Optimisation of a Pressurised Liquid Extraction Method for Haloanisoles in Cork Stoppers. *Anal. Chim. Acta.* **2005**, *540*, 17-24.
- Draisci, R.; Marchiafava, C.; Palleschi, L.; Cammarata, P.; Cavalli, S. Accelerated Solvent Extraction and Liquid Chromatography-Tandem Mass Spectrometry Quantitation of Corticosteroid Residues in Bovine Liver. *J. Chromatog., B.: Anal. Technol. Biomed. Life Sci.* **2001**, *753*, 217-223.
- Gentili, A.; Perret, D.; Marchese, S.; Sergi, M.; Olmi, C.; Curini, R. Accelerated Solvent Extraction of Confirmatory Analysis of Sulfonamide Residues in Raw Meat and Infant Foods by Liquid Chromatography Electrospray Tandem Mass Spectrometry. *J. Agric. Food Chem.* **2004**, *52*, 4614-4624.
- Dionex Corporation (now part of Thermo Fisher Scientific). Extraction and Clean Up of Acrylamide in Complex Matrices Using ASE Followed by Liquid Chromatography, Tandem Mass Spectrometry (LC-MS/MS). Application Note 358; LPN 1871. Sunnyvale, CA, 2006.
- Poerschmann, J.; Carlson, R. New Fractionation Scheme for Lipid Classes Based on "In-Cell Fractionation" Using Sequential Pressurized Liquid Extraction. *J. Chromatog., A.* **2006**, *1127*, 18-25.
- Poerschmann, J.; Trommler, U.; Biedermann, W.; Truyen, U.; Lucker, E. Sequential Pressurized Liquid Extraction to Determine Brain-Originating Fatty Acids in Meat Products as Markers in Bovine Spongiform Encephalopathy Risk Assessment Studies. *J. Chromatog., A.* **2006**, *1127*, 26-33.

www.thermoscientific.com/dionex

©2012 ThermoFisher Scientific Inc. All rights reserved. ISO is a trademark of the International Standards Organization. Florisil is a registered trademark of U.S. Silica Co. Isolute is a registered trademark of Biotage AB. All other trademarks are the property of ThermoFisher Scientific Inc. and its subsidiaries. This information is presented as an example of the capabilities of ThermoFisher Scientific Inc. products. It is not intended to encourage use of these products in any manner that might infringe the intellectual property rights of others. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.



Thermo Scientific Dionex products are designed, developed, and manufactured under an ISO 9001 Quality System.

Australia +61 3 9757 4486	China +852 2428 3282	Ireland +353 1 644 0064	Sweden +46 8 473 3380
Austria +43 1 616 51 25	Denmark +45 36 36 90 90	Italy +39 02 51 62 1267	Switzerland +41 62 205 9966
Benelux +31 20 683 9768	France +33 1 39 30 01 10	Japan +81 6 6885 1213	Taiwan +886 2 8751 6655
+32 3 353 42 94	Germany +49 6126 991 0	Korea +82 2 3420 8600	UK +44 1276 691722
Brazil +55 11 3731 5140	India +91 22 2764 2735	Singapore +65 6289 1190	USA and Canada +847 295 7500

Thermo
SCIENTIFIC

Part of Thermo Fisher Scientific